




This accessory is intended to provide mobility and assist in transferring liquid nitrogen from the model LD25. The easy-to-assemble stand has a 5 wheel cast-aluminum dolly for handling convenience and side brackets to support the weight of the Dewar while you tip the unit forward to pour. When not in use, the weight of the contents returns the Dewar to an upright position. To position the Dewar in the stand, secure the unit in the pivot ring so that it balances itself upright when in storage or travel and yet allows you to safely control the unit while pouring.

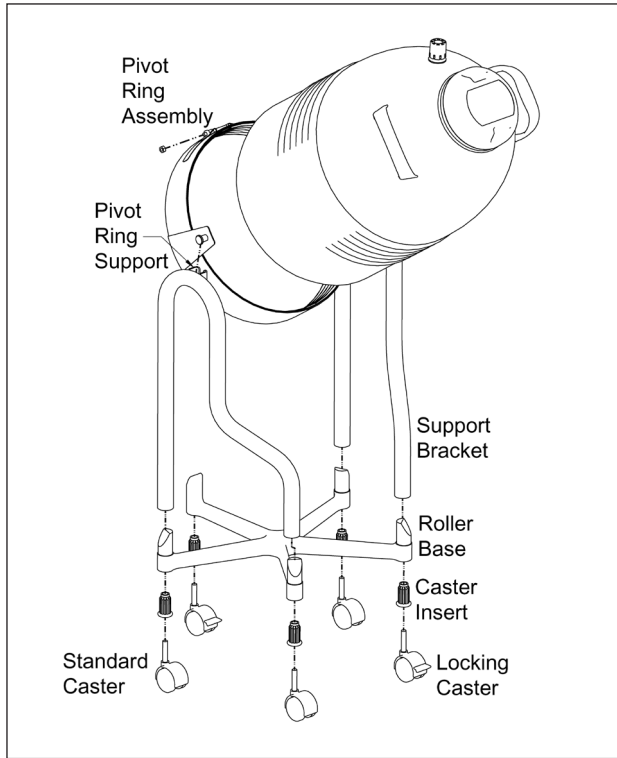

Tipping Stand Drawing

Caution:

Avoid liquid nitrogen over-filling. Adhesives that are subjected to extreme cold should be inspected and, if needed, should be replaced.

LIQUID NITROGEN CAN PRESENT HAZARDS THAT RESULT IN PERSONAL INJURY. To protect yourself and others, read "Handle with Care", and the Operating Instructions that are provided with Dewar.

Assembly Instructions

- Slide the support brackets down, over the smooth roller base pegs, so that the pivot ring supports are directed inward.
- Remove the nut on the pivot ring to loosen the assembly. Position the Dewar in place, leaving one corrugation showing above the ring. Tighten the nut back into place.
- Slide the pivot ring assembly into place and press down into position. If necessary, reposition the Dewar in the pivot ring to get the proper balance described above.

REPLACEMENT PARTS

| Description | Part Number |
|------------------------|-------------|
| Roller Base | 366785 |
| Support Bracket, Left | 366584 |
| Support Bracket, Right | 366585 |
| Caster, Standard | 367567 |
| Caster, Locking | 367568 |
| Pivot Ring Assembly | 366619 |
| Caster Insert | 367306 |
| Rubber Grip, Contoured | 366583 |

CONTACT US

US Toll Free: 844-273-7517

Global: +43 7485 606 286 / +1 614-438-7968

Sales Support/Order Placement:
customerservice@worthingtonindustries.com

Technical Services:
techservices@worthingtonindustries.com

